

Sample Preparation Kit (96 reactions)

PROTOCOL – Pelleted cells & precipitated protein



INTRODUCTION

Sample preparation is one of the essential steps of bottom-up proteomics. The PreOmics sample preparation kit is designed to assist you achieving best results with few sample preparation steps and little hands-on time. For sample-specific protocols and optimization contact us or visit our website at www.preomics.com.

KIT CONTENTS

The kit contains all you need to perform a complete sample preparation. It includes all chemicals to denature, reduce and alkylate proteins as well as the enzymes to perform a tryptic digest and a final clean up.

COMPONENT	CAP	QUANTITY	BUFFER PROPERTIES				DESCRIPTION	STORAGE
			Organic	Acidic	Basic	Volatile		
DIGEST	red	24x					Enzyme <i>Trypsin</i> -mix to digest proteins.	-20 °C
RESUSPEND	yellow	4x 2 mL				✓	Protease reconstitution buffer for enzymes.	RT
LYSE	brown	12x 1 mL				✓	Denature, reduce and alkylate proteins.	RT
STOP	black	12x 1 mL	✓	✓		✓	Stop the enzymatic activity.	RT
WASH 1	blue	12x 2 mL	✓	✓		✓	Clean up peptides from hydrophobic contaminants.	RT
WASH 2	green	12x 2 mL				✓	Clean up peptides from hydrophilic contaminants.	RT
ELUTE	violet	12x 2 mL	✓			✓	Elute the peptides from the cartridge.	RT
LC-LOAD	white	12x 1 mL				✓	Load peptides on reversed-phase LC-MS column.	RT
CARTRIDGES		96x					Cartridges for 1 to 100 µg protein starting material.	RT
WASTE PLATE		1x					Deep well plate for collecting waste after washes.	RT
MTP PLATE		1x					LoBind plate for collecting peptides after elution.	RT
ADAPTER PLATE		1x					Enables placing cartridges in the 96well MTP plate.	RT
CAPS		96x					Caps to optionally close the cartridge's bottom.	RT


PRE-REQUISITES

Common lab equipment is required for the sample preparation.

EQUIPMENT	QUANTITY AND DESCRIPTION
PIPETTE	Careful sample handling and pipetting reduces contaminations and improves quantification.
SAMPLE	Pelleted cells or precipitated protein. For other materials ask us for adapted protocols.
96x-WELL	96 deep well and PCR plates with skirt to balance WASTE & MTP plates in centrifuge.
HEATING BLOCK*	2x MTP plate heaters are recommended to help protein denaturation and during digestion.
CENTRIFUGE	Wing-bucket MTP plate centrifuges are necessary for loading, washing and elution.
SONICATOR	If you have DNA in your sample, shear it by sonication (e.g. Diagenode Bioruptor®).
SPEED-VAC	Vacuum manifolds evaporate volatile buffers from the eluate before LC-MS.

* a heating shaker adapter for MTP plate shakers can be purchased from PreOmics

PROCEDURE

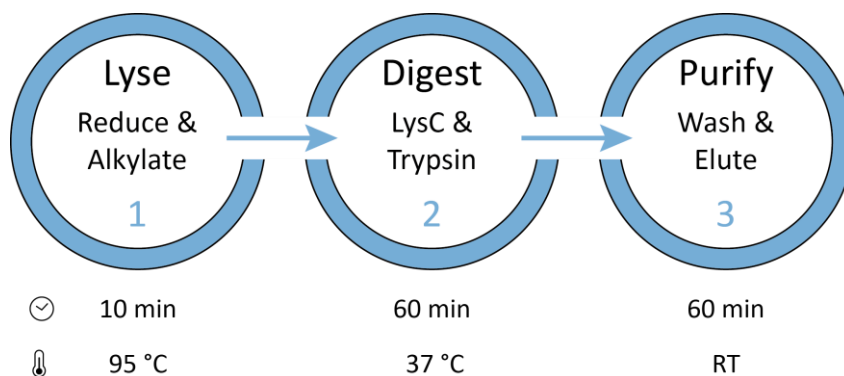

Sample Preparation Kit (96 reactions)

PROTOCOL – Pelleted cells & precipitated protein



PROTOCOL

1. LYSE

- 1.1. Add 50 µL **LYSE** ● to 1 - 100 µg of sample and place it in a pre-heated HEATING BLOCK (95 °C; 1,000 rpm; 10 min).
- 1.2. Optional: Spin down droplets (RT; max. 300 rcf; 10 sec).
- 1.3. If you expect DNA, shear sample in a SONICATOR (10 cycles 30 sec ON/OFF).
- 1.4. Transfer sample to **CARTRIDGE** and cool down (RT). Be careful not to damage the bottom layer of **CARTRIDGE**.
- 1.5. Use **ADAPTER PLATE** to place **CARTRIDGES** on top of the **WASTE PLATE** tube. Label all plate and wells.

2. DIGEST

- 2.1. Add 210 µL **RESUSPEND** ● to **DIGEST** ● (1 tube for 4 reactions), shake (RT; 500 rpm; 10 min), pipette up/down.
- 2.2. Add 50 µL **DIGEST** ● to **CARTRIDGE** and place it in a pre-heated HEATING BLOCK (37 °C; 500 rpm; 1-3 hours).
- 2.3. Add 100 µL **STOP** ● to **CARTRIDGE** (precipitation may happen), shake (RT; 500 rpm; 1 min), pipette up/down. *SP*

3. PURIFY

- 3.1. Spin **CARTRIDGE** in a CENTRIFUGE (3,800 rcf; 1-3 min). If needed, adjust values to ensure complete flow-through.
- 3.2. Add 200 µL **WASH 1** ● to **CARTRIDGE**, repeat step 3.1., discard flow-through.
- 3.3. Add 200 µL **WASH 2** ● to **CARTRIDGE**, repeat step 3.1., discard flow-through. *SP*
- 3.4. Use **ADAPTER PLATE** to place **CARTRIDGES** on top of the **MTP PLATE**. Label plate and wells.
- 3.5. Add 100 µL **ELUTE** ● to **CARTRIDGE**, repeat step 3.1., keep flow-through in **MTP PLATE**.
- 3.6. Repeat step 3.5., keep flow-through in the same **MTP PLATE**.
- 3.7. Remove **CARTRIDGES** and place **MTP PLATE** in a SPEED-VAC (45 °C; until completely dry).
- 3.8. Add **LC-LOAD** ○ to **MTP PLATE**. Aim for 1 g/L concentration (e.g. 100 µL to 100 µg protein starting material).
- 3.9. Sonicate **MTP PLATE** in a SONICATOR (10 cycles 30 sec ON/OFF). *SP*

SP - Storage Point: At this point, close the peptide containing tube or **CARTRIDGE** (use a **CAP** for bottom). Peptides can be frozen at -20 °C. Storage of peptides should not exceed 2 weeks at -20 °C. For extended storage, finish the protocol and store at -80 °C.

DATA ANALYSIS

Consider the following as fixed modifications in your database search:

MODIFICATION	DESCRIPTION	COMPOSITION	SPECIFICITY	MASS	UNIMOD #
ALKYLATION	Carbamidomethyl on Cysteine	C ₂ H ₃ NO	[C]	+57 Da	4

DISCLAIMER

Products are warranted to operate or perform substantially in conformance with published Product specifications in effect at the time of sale, as set forth in the Product documentation, specifications and/or accompanying package inserts ("Documentation"). No claim of suitability for use in applications regulated by FDA is made. The warranty provided herein is valid only when used by properly trained individuals. Unless otherwise stated in the Documentation, this warranty is limited to three months from date of shipment when the Product is subjected to normal, proper and intended usage. This warranty does not extend to anyone other than Buyer. Any model or sample furnished to Buyer is merely illustrative of the general type and quality of goods and does not represent that any Product will conform to such model or sample. NO OTHER WARRANTIES, EXPRESS OR IMPLIED, ARE GRANTED, INCLUDING WITHOUT LIMITATION, IMPLIED WARRANTIES OF MERCHANTABILITY, FITNESS FOR ANY PARTICULAR PURPOSE, OR NON-INFRINGEMENT. BUYER'S EXCLUSIVE REMEDY FOR NONCONFORMING PRODUCTS DURING THE WARRANTY PERIOD IS LIMITED TO REPAIR, REPLACEMENT OF OR REFUND FOR THE NONCONFORMING PRODUCT(S) AT SELLER'S SOLE OPTION. THERE IS NO OBLIGATION TO REPAIR, REPLACE OR REFUND FOR PRODUCTS AS THE RESULT OF (I) ACCIDENT, DISASTER OR EVENT OF FORCE MAJEURE, (II) MISUSE, FAULT OR NEGLIGENCE OF OR BY BUYER, (III) USE OF THE PRODUCTS IN A MANNER FOR WHICH THEY WERE NOT DESIGNED, OR (IV) IMPROPER STORAGE AND HANDLING OF THE PRODUCTS. Unless otherwise expressly stated on the Product or in the documentation accompanying the Product, the Product is intended for research only and is not to be used for any other purpose, including without limitation, unauthorized commercial uses, in vitro diagnostic uses, ex vivo or in vivo therapeutic uses, or any type of consumption by or application to humans or animals.